

Vitamin Analyzer

Vitamins are unstable compounds, which are easily oxidized and destroyed during sample preparation procedures. For that reason, a simple assay method that reduces degradation of vitamins needs to be established in order to produce accurate results.

Typically, vitamins have been analyzed by spectrophotometric and microbiological procedures, which involve lengthy processes due to complicated sample preparation procedures, and often result in poor reproducibility.

Recently, HPLC has been used for analyses of vitamins in order to overcome these problems, and produce accurate and reproducible results.

YL Vitamin Analyzer is dedicated to analyze water-soluble and fat-soluble vitamins.

• Useful Information

• Water-soluble vitamin

Vitamin C (Ascorbic acid), B1 (Thiamine), B2 (Riboflavin), B3 (Niacin), B6 (Pyridoxine), B12 (Cyanocobalamin), Folic acid, Pantothenic acid


• Fat-soluble vitamin

Vitamin A (Retinol), D, E, K (K1, K2, K3)


Sample preparation procedures and the analysis of Vitamins

(1) Sample preparation procedure of Vitamins

• Sample preparation procedure for water-soluble vitamins

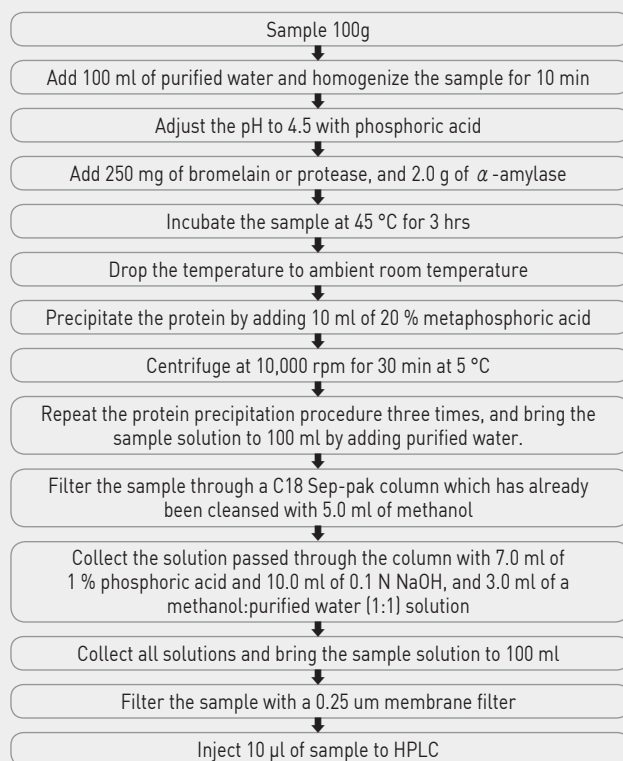

• Sample preparation procedure for fat-soluble vitamins

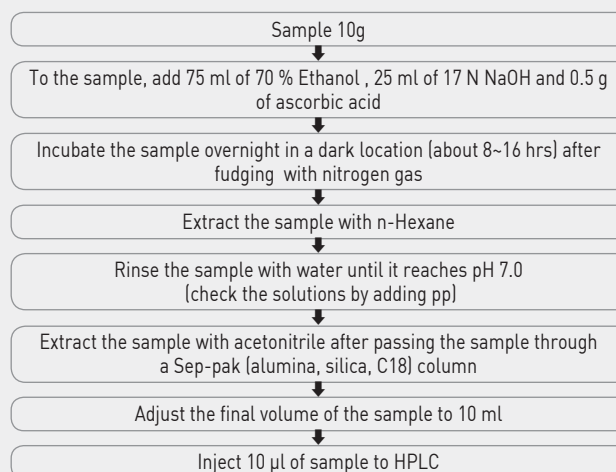

• Recommended Columns

- Water-soluble vitamins : Aqueous C8 (4.6 mm, 150 mm)
- Fat-soluble vitamins : C8 (4.6 mm, 150 mm)

• Application

- Vitamins in powdered milk
- Vitamins in beverage
- Vitamins in natural foods
- Vitamins in a medicine promoting nutrition

Алматы (7273)495-231
 Ангарск (3955)60-70-56
 Архангельск (8182)63-90-72
 Астрахань (8512)99-46-04
 Барнаул (3852)73-04-60
 Белгород (4722)40-23-64
 Благовещенск (4162)22-76-07
 Брянск (4832)59-03-52
 Владивосток (423)249-28-31
 Владикавказ (8672)28-90-48
 Владимир (4922)49-43-18
 Волгоград (844)278-03-48
 Вологда (8172)26-41-59
 Воронеж (473)204-51-73
 Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
 Ижевск (3412)26-03-58
 Иркутск (395)279-98-46
 Казань (843)206-01-48
 Калининград (4012)72-03-81
 Калуга (4842)92-23-67
 Кемерово (3842)65-04-62
 Киров (8332)68-02-04
 Коломна (4966)23-41-49
 Кострома (4942)77-07-48
 Краснодар (861)203-40-90
 Красноярск (391)204-63-61
 Курск (4712)77-13-04
 Курган (3522)50-90-47
 Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
 Москва (495)268-04-70
 Мурманск (8152)59-64-93
 Набережные Челны (8552)20-53-41
 Нижний Новгород (831)429-08-12
 Новокузнецк (3843)20-46-81
 Ноябрьск (3496)41-32-12
 Новосибирск (383)227-86-73
 Омск (3812)21-46-40
 Орел (4862)44-53-42
 Оренбург (3532)37-68-04
 Пенза (8412)22-31-16
 Петрозаводск (8142)55-98-37
 Псков (8112)59-10-37
 Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
 Рязань (4912)46-61-64
 Самара (846)206-03-16
 Санкт-Петербург (812)309-46-40
 Саратов (845)249-38-78
 Севастополь (8692)22-31-93
 Саранск (8342)22-96-24
 Симферополь (3652)67-13-56
 Смоленск (4812)29-41-54
 Сочи (862)225-72-31
 Ставрополь (8652)20-65-13
 Сургут (3462)77-98-35
 Сыктывкар (8212)25-95-17
 Тамбов (4752)50-40-97
 Тверь (4822)63-31-35

Тольятти (8482)63-91-07
 Томск (3822)98-41-53
 Тула (4872)33-79-87
 Тюмень (3452)66-21-18
 Ульяновск (8422)24-23-59
 Улан-Удэ (3012)59-97-51
 Уфа (347)229-48-12
 Хабаровск (4212)92-98-04
 Чебоксары (8352)28-53-07
 Челябинск (351)202-03-61
 Череповец (8202)49-02-64
 Чита (3022)38-34-83
 Якутск (4112)23-90-97
 Ярославль (4852)69-52-93


Water Soluble Vitamins

- Mobile phase : A- Ultrapure water (10 mM NaH₂PO₄ + 10 mM Na₂HPO₄ + H₃PO₄ 0.1 %)
B- MeOH

Gradient program


Time (min)	Flow Rate (ml/min)	Mobile Phase	
		A (%)	B (%)
0	1.0	100	0
2	1.0	100	0
15	1.0	40	60
16	1.0	40	60

- Column oven : 35 °C
- Detector : UV/Vis 280 nm
- Column : Atlantis dC18 (4.6 mm, 150 mm, 5 µm)
- Injection volume : 20 µl sample loop

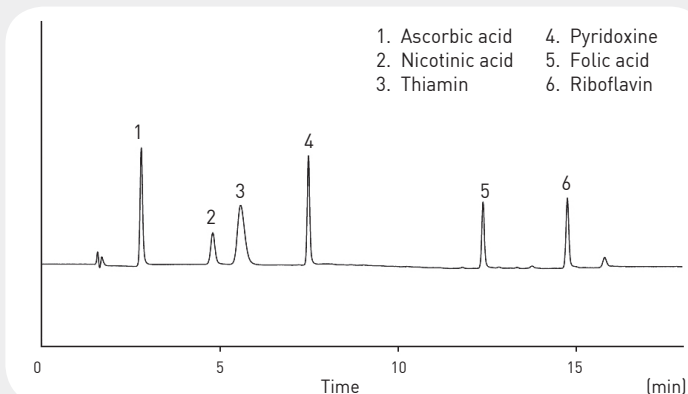


Fat Soluble Vitamins

- Mobile phase : ACN:MeOH = 75:25
- Flow rate : 1.0 ml/min
- Column oven : 35 °C
- Detector : UV/Vis 280 nm
- Column : C18 (4.6 mm, 250 mm, 5 µm)
- Injection volume : 20 µl sample loop

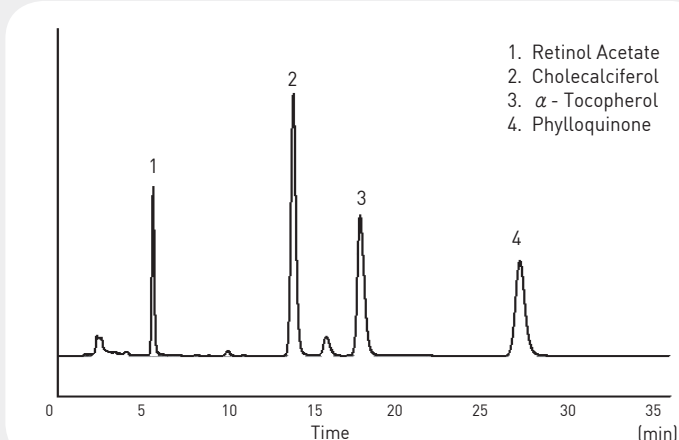

Fat Soluble Vitamins E (α - Tocopherol, Tocopherol Acetate)

- Mobile phase : MeOH 49 % : Water 1 %
- Flow rate : 1 ml
- Column oven : 25 °C
- Detector : UV 284 nm
- Column : Xbridge C18 (4.6 mm, 150 mm, 5 µm)

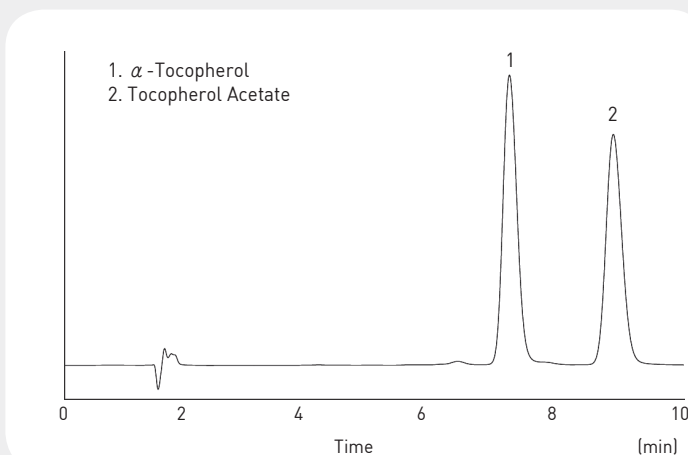

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Саранск (8342)22-96-24
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93